


Fold back and forth several times along joint lines before assembling.

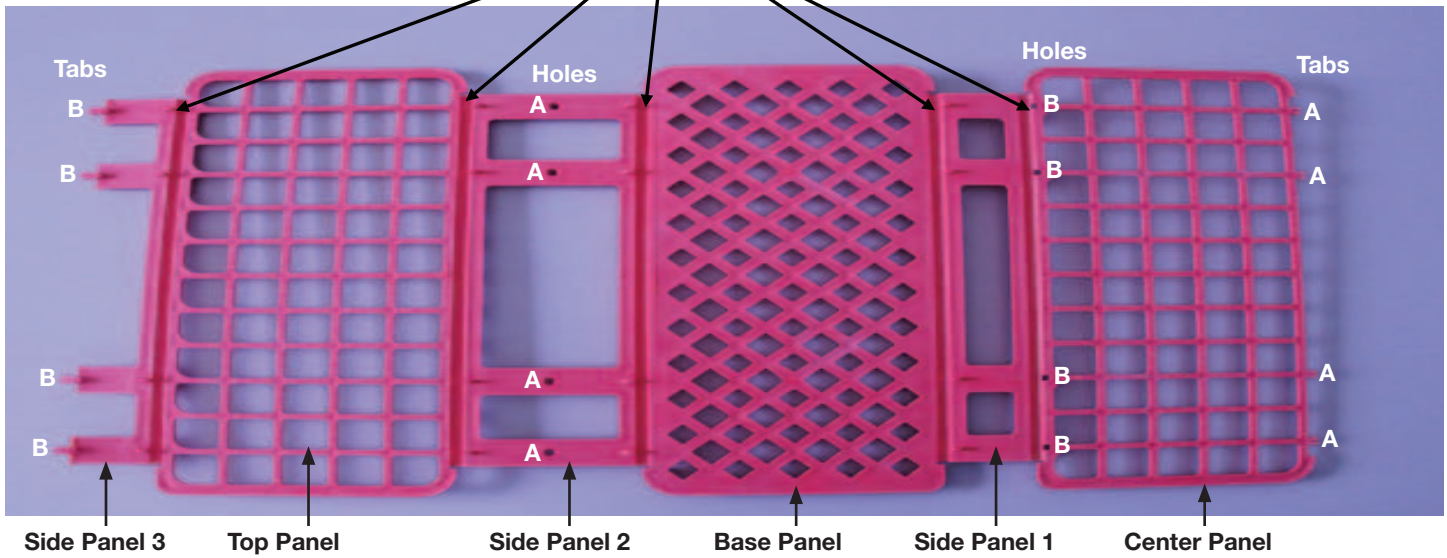

To Assemble:

1. Fold rack back and forth several times along joint lines before assembling.
2. Fold Center Panel up and over until parallel with Base Panel.
3. Fold Side Panel 2 up and align Tabs A with Holes A. Snap into place. **Note:** It is easiest to stand rack sideways on table or workbench and snap tabs into holes one at a time using equal pressure on left and right side of each tab.
4. Fold Top Panel over until parallel with Base and Center Panels.
5. Lay the partially assembled rack flat on table or workbench. Fold Side Panel 3 down and snap Tabs B into Holes B along edge of Center Panel. Again, snap tabs one at a time using equal pressure on both sides of the tabs simultaneously.

